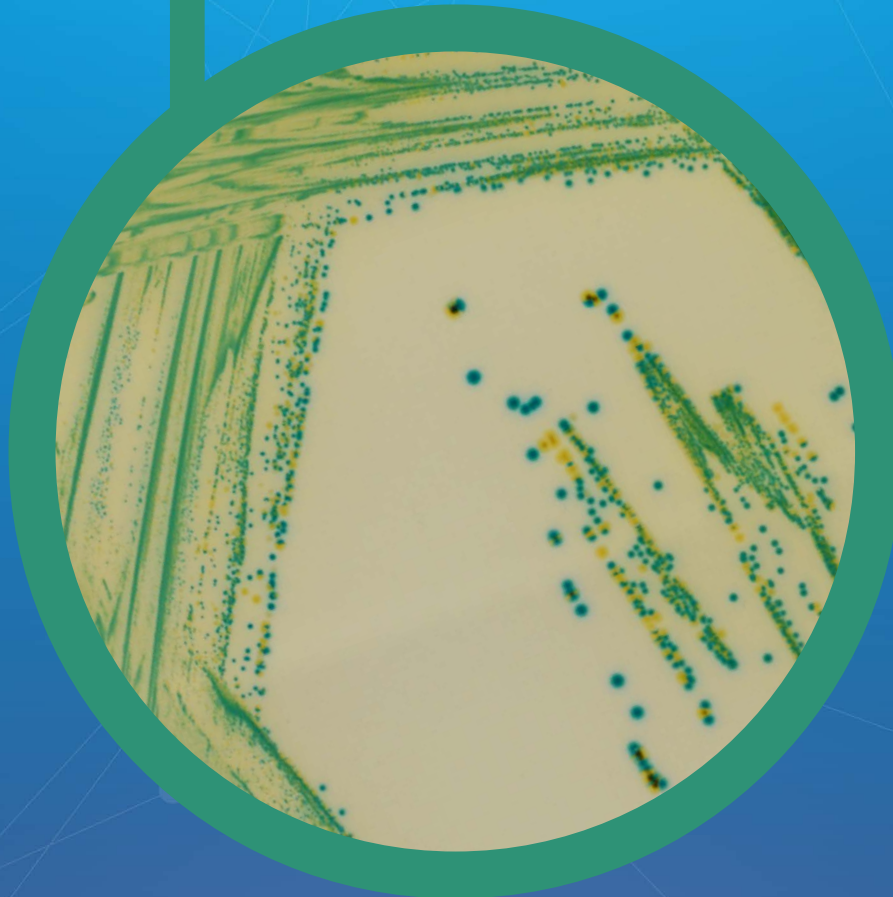
# Leaflet

LF-EXT-065

Version 4.0

## ● CHROMagar™ Cronobacter

Click below:

EN

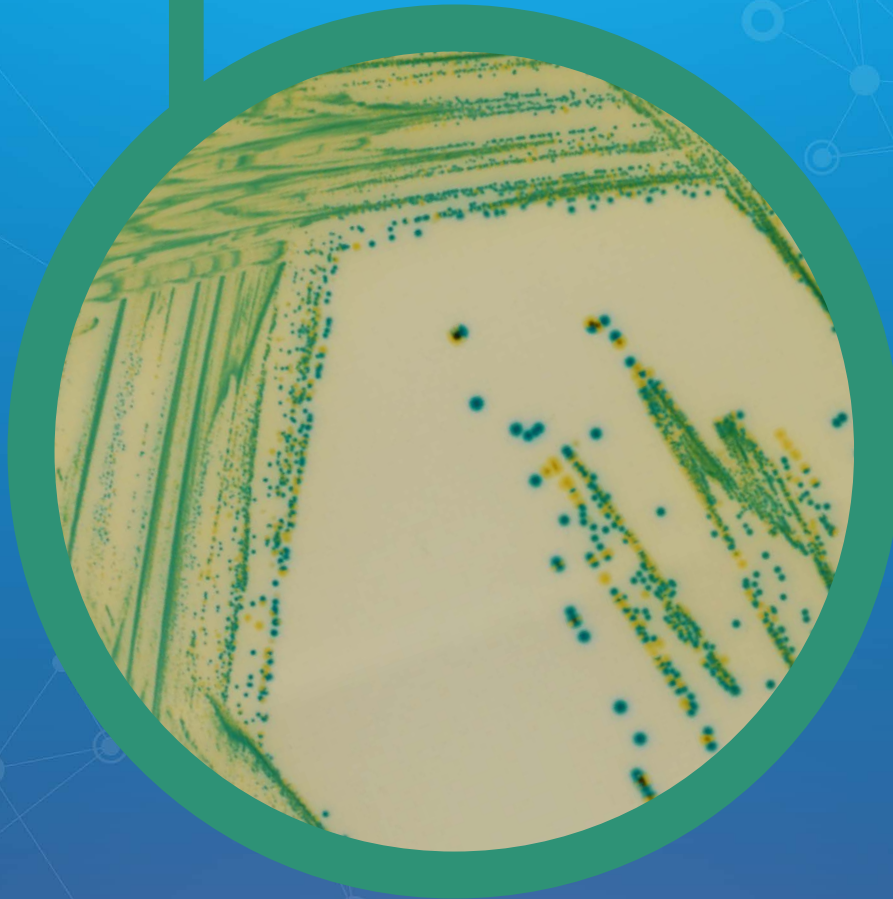
FR

ES


CHR  Magar™  
The Chromogenic Media Pioneer

**Water industry**

# ● CHROMagar™ Cronobacter

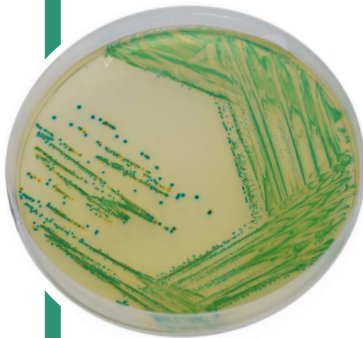**Chromogenic Cronobacter Isolation (CCI) agar,  
in accordance with ISO 22964 standard**

## Plate Reading

- *Cronobacter* spp.  
→ green to blue
- Other Gram (-)  
→ colourless, clear green, black, yellow
- Gram (+) bacteria  
→ inhibited

## Chromogenic Cronobacter Isolation (CCI) agar, in accordance with ISO 22964 standard

### Background

«*Cronobacter* is an ubiquitous telluric microorganism found in water, soil, plants, dust and many other living things. *Cronobacter* has been isolated from many foods of plant or animal origin, both dehydrated, smoked, frozen, fermented, raw and cooked.»

Extracts from ANSES, April 2017

CHROMagar™ Cronobacter is the Chromogenic Cronobacter Isolation (CCI) agar, manufactured in accordance with ISO 22964 standard.

### Medium Performance

**1 ISO 22964 STANDARD**  
CHROMagar™ Cronobacter is a chromogenic medium for detection of *Cronobacter* spp. in food, animal feed and environmental samples according to the ISO 22964 standard.

**2 INTENSE GREEN TO BLUE COLONIES**  
*Cronobacter* spp. will grow in green to blue colonies while most other bacteria will be inhibited or grow in colourless colonies, after Incubation at 41.5 °C for 24 h.

### Medium Description

Powder Base	
Total .....	32.7 g/L
Agar .....	15.3
Casein Peptone .....	7.0
Yeast extract .....	3.0
Sodium chloride .....	5.0
5-bromo-4-chloro-3-indolyl- $\alpha$ -D-glucopyranoside ...	0.15
Sodium deoxycholate .....	0.25
Ammonium iron(III) citrate .....	1.0
Sodium thiosulfate .....	1.0
Storage at 15/30 °C - pH: 7.3 +/- 0.2	
Shelf Life .....	2 years

Usual Samples	Food, animal feed, environmental samples.
Procedure	Appropriate pre-enrichment and enrichment steps. Streaking 10 $\mu$ L. Incubation 24 h at 41.5 °C Aerobic conditions.

Please read carefully the instructions for use (IFU document) available on [www.CHROMagar.com](http://www.CHROMagar.com)

## Order References

Please use these product references when contacting your local distributor:

500 G pack ..... MC123-500G

Manufacturer: CHROMagar, 29 avenue George Sand,  
93210 La Plaine Saint-Denis - France

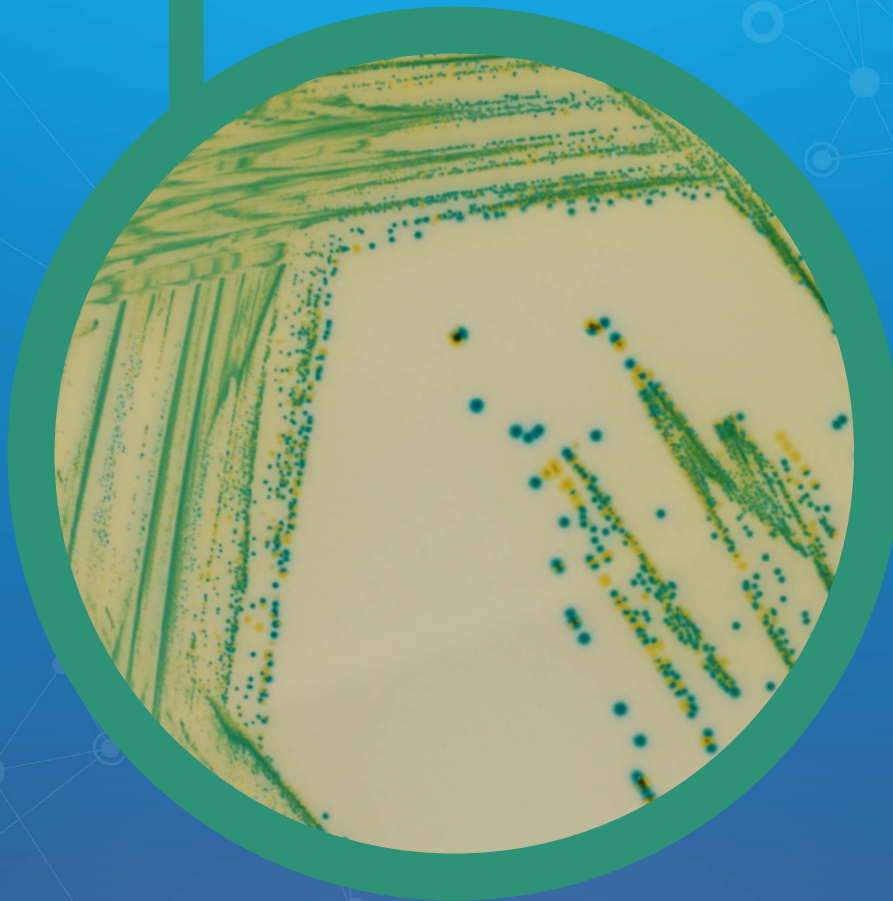
Email: [CHROMagar@CHROMagar.com](mailto:CHROMagar@CHROMagar.com)

Website: [www.CHROMagar.com](http://www.CHROMagar.com)

Find your nearest distributor on  
[www.CHROMagar.com/contact](http://www.CHROMagar.com/contact)

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